

PDQ™ Protease Assay

Catalog Numbers: 0201

Product Description

The **PDQ™ (Protease Determine Quick) Protease Assay** is a colorimetric assay used to detect protease activity in aqueous samples. The proprietary substrate responds to a wide range of proteases including serine, metallo, aspartate and cysteine proteases such as collagenase, proteinase K, papain, pepsin, bromelin, ficin, trypsin and chymotrypsin. PDQ™ can be used with just a few simple steps to measure protease activity and requires no centrifugation. The substrate is a cross-linked matrix containing protein substrate and a dye-protein conjugate. Protease activity is detected spectrophotometrically with increasing optical density proportional to increasing enzyme activity and can detect nanogram quantities.

Instructions for Use

1. Allow the vials to come to room temperature before running the assay. Pour off storage solution.
2. To construct a standard curve, prepare a ten-fold dilution series of the control protease. This control solution should contain the protease for which the test is being run. Trypsin is supplied as a general protease control. The trypsin can be used in a ten-fold dilution series; e.g., 400 to 0.4 BAEE units/ml (280 - 0.28 µg/ml). Any buffer can be used with the PDQ™ matrix. (See Technical Considerations). The standard trypsin reaction buffer is 10mM Tris-Cl, pH 8.0.
3. Add 0.5 ml of each control solution to duplicate vials.
4. Prepare test samples containing putative protease activity in the same buffer as the control protease. Add 0.5 ml of the test sample to duplicate vials.
5. Replace caps onto vials.
6. Incubate vials at 37°C for 1 h. (A shorter or longer incubation time may be required. See Technical Considerations).
7. Add 0.5 ml 0.2N NaOH to each vial to stop the reaction and amplify color. Either gently swirl the vial or replace the cap and invert the vial once to mix contents.
8. Transfer the contents (being careful not to disturb the semi-solid matrix) to a standard 1 ml cuvette. (Transfer 0.2 ml of liquid if using a microplate reader.)
9. Record the absorbance at 450 nm.
10. Average the results from duplicate vials and use the control samples to generate a standard curve, plotting concentration (or amount) of control protease vs. A450. Compare test sample results against the standard curve to obtain the amount of protease activity.

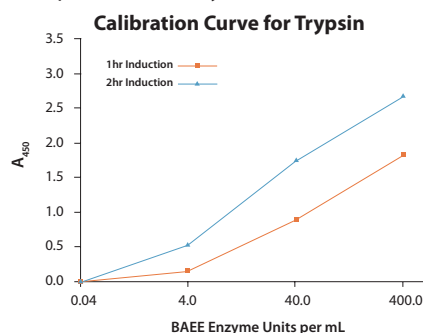


Figure. Calibrator curves generated using the PDQ™ Colorimetric Assay. Trypsin was at 400, 40, 4.0, and 0.4 BAEE Units/mL incubated for 1 or 2 hours.



Athena Enzyme Systems™
1450 South Rolling Road
Baltimore, MD 21227
USA
AthenaES®
a division of Athena Environmental Sciences, Inc.

T (MD): 410-455-6319
T (USA): 888-892-8408
F: 410-455-1155
aesinfo@athenaes.com

Technical Considerations

1. Buffers that can be used to prepare test samples may range in pH from 4 to 10.
2. The amount of time required for incubation can be varied to optimize the assay. Up to 24 hr incubations are possible which will increase the detection limits of the assay. Alternatively, for applications with high amounts of enzyme, 15 min may suffice. The exact conditions should be determined empirically.
3. As with any assay, it is necessary to run a standard curve with control protease every time the assay is performed with unknown protease samples.
4. When using the vials, the stopped reaction contents must be transferred to a cuvette to measure the absorbance. The transfer of the contents away from the PDQ™ matrix should be done as soon as possible after the addition of NaOH, since the NaOH will degrade the matrix over time. However, the contents can be left on the matrix for up to one hour before transfer without compromising the accuracy of the test.
5. Once the stopped reaction contents are transferred away from matrix, they can be left for up to 24 hrs covered (to keep out light) at room temperature before measuring the absorbance without loss of signal.
6. To use the Colorimetric Assay for opaque samples, run the reaction as described. Remove the reaction liquid and then back digest the remaining substrate matrix with Trypsin, Subtilisin or Proteanase K. The signal generated from the back digestion is inversely proportional to the amount of enzyme activity.

Product Specifications

0201 Colorimetric PDQ™ Protease Assay	
Kit Size	48 vials of 200µL of pre-made matrix
Control	Trypsin supplied as control: 0.5mL of 0.7mg/mL (1420 BAEE units/mg) in reaction buffer (10mM Tris-Cl, pH 8.0).
Storage	4°C
Stability	3 months

Materials Required but not Provided

0.2N NaOH (for colorimetric assay only)
Spectrophotometer for reading absorbance at 450nm (standard assays)